DS1070
055006

EN	FR
IT	EL
BG	SL
RO	HR

v2.2

WWW.NIKOLAOUTOOLS.COM

GENERAL INSTRUCTIONS

Attention: Read the manual carefully before use. Follow all the instructions contained in the manual in order to keep the scale in good working order. Save the manual in a safe place for future reference.

- Never place loads on the scale that exceed its weighing capacity. Overloading the scale will damage it.
- Avoid exposing the unit to extreme heat or cold. The scale works best when operated at normal room temperature.
- Always allow the scale to acclimate to normal room temperature for at least 1 hour before use.
- Give the scale sufficient warm up time. In most cases, 30 seconds of warm up time suffice to ensure that the internal components have stabilised before calibration.
- Keep the area in which the scale is placed clean. Dust, dirt, moisture, vibration, air currents and a close distance to other electronic equipment can all cause an adverse effect on the reliability and accuracy of the scale.
- Handle the scale with care. Gently place all items to be weighed onto the platform. Do not abuse the scale as this may permanently damage the internal sensor.
- Avoid shaking, dropping or hitting the scale. This is a precision instrument and must be handled with extreme care.
- Only operate the scale on a stable surface, free of vibration.

OPERATING INSTRUCTIONS

Basic operation

1. Place the scale on a horizontal, flat surface. Press the [ON/OFF] button to turn on the scale.
2. Wait until "0.0" is displayed.
3. Put the item(s) on the weighing platform.
4. Using the [MODE] button, you can switch between the weighing units supported by the scale.
5. Using the [LIGHT] button, you can turn on or off the back light.

Tare function

1. Turn on the scale as described above.
2. Place an empty container on the platform.
3. Press the [TARE] button and wait until "0.0" is displayed.
4. Add items into the empty container.

Counting function

1. Turn on the scale and wait until "0.0" is displayed on the screen.
2. Place a sample amount (25pcs, 50pcs, 75pcs or 100pcs) of the item you wish to weigh and count onto the platform.
3. Press and HOLD the [PCS] button for 2 seconds. The screen will display a random number, for example "100".
4. Press the [MODE] button to select the sample amount of the item that you have put onto the platform.
5. Press the [PCS] button to confirm the above information. The screen will now display "pcs" in the corner, indicating that the scale is in counting mode.
6. Remove the sample quantity from the platform.
7. You now can place any amount of item(s) on the scale and the screen will display a count.

Note: To switch the scale to weighing mode, press the [MODE] button.

Note: To switch the scale to counting mode, press the [PCS] button.

Calibration

1. Reset the scale by pressing the [ON/OFF] button.
2. Wait until "0.0" is displayed on the screen and then press and HOLD the [MODE] button for 3 seconds. Once the "CAL" indication is displayed on the screen, release the [MODE] button.
3. Press the [MODE] button again. The "CAL" indication will be displayed on the screen followed by the required calibration weight.
4. Add the required calibration weight. The "PASS" indication will be displayed on the screen and the scale will return to normal weighing mode.
5. Remove the required calibration weight and press the [ON/OFF] button to turn off the scale. Calibration is now finished.

TECHNICAL SPECIFICATIONS

Technical Data

Technical Data	
Model	DS1070
Weighing capacity	2.000 gr
Resolution	0.1 gr

* This product is intended for laboratory use only.

* The manufacturer reserves the right to make minor changes to product design and technical specifications without prior notice unless these changes significantly affect the performance and safety of the products. The parts described / illustrated in the pages of the manual that you hold in your hands may also concern other models of the manufacturer's product line with similar features and may not be included in the product you just acquired.

* To ensure the safety and reliability of the product and the warranty validity, all repair, inspection, repair or replacement work, including maintenance and special adjustments, must only be carried out by technicians of the authorized service department of the manufacturer.

* Always use the product with the supplied equipment. Operation of the product with non-provided equipment may cause malfunctions or even serious injury or death. The manufacturer and the importer shall not be liable for injuries and damages resulting from the use of non-conforming equipment.

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